Méthode ESAO[®]

ESAO[®] method

**Ref :
453 011**

Français – p 1

English – p 3

Version : 6006

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1 General information

Manufactured like an Initio type case, the synchro trigger system helps synchronise a measurement during a mechanical action (with ESAO®4, ESAO®3.1 or VTT interface).

It helps, for example, synchronise the measurement during the simulation of a seismic wave with VTT and the simplified Ex.A.O. seismograph (**Part no. 453 010 79**).

It is sturdy and can bear an impact of 5 joules, i.e. the fall of a 500 g ball from a height of 1 metre.

2 Use with ESAO®

- With 2 cables equipped with safety sockets, connect the trigger system to the ESAO® console using the Synchro input (red and black sockets).
- Set up the software to synchronise the measurement on the high-to-low transition synchro channel.
- Start the acquisition.
- Using a mallet or any other object that enables the application of an impact of moderate intensity (less than 5 joules), hit on the central metal portion of the trigger system.

3 Use with VTT®

The synchro trigger system may be used with VTT.

**In this case, the VTT console must be equipped with version 1.15 or later of the recording software.
You can find an updated version of this software available for free download on the Internet or you may ask us for it by calling our technical support.**

- With 2 cables equipped with safety sockets, connect the trigger system to the VTT console using the Synchro input (red and black sockets on the side of the VTT console).
- Set up the recording software to synchronise the measurement.
- Start the acquisition.
- Using a mallet or any other object that enables the application of an impact of moderate intensity (less than 5 joules), hit on the central metal portion of the trigger system.

4 Technical specifications

- Contact by metallic grounding
- 2 outputs per safety plug
- Maximum operating voltage 12V/200mA.

5 After-sales service

The device is under a 2-year guarantee, it must be sent back to our workshops.

For any repairs, adjustments or spare parts, please contact:

JEULIN – TECHNICAL SUPPORT
Rue Jacques Monod
BP 1900
27 019 EVREUX CEDEX FRANCE
+33 (0)2 32 29 40 50

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SA capital 3 233 762 € - Siren R.C.S. B 387 901 044 - Siret 387 901 04400017

Direct connection for technical support

A team of experts at your disposal from Monday to Friday (opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediatly to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

Free service * :
+ 33 (0)2 32 29 40 50

** Call cost not included*