Tooxy

Capteur

Tooxy

Sensor

Ref :
482 201

Français – p 1

English – p 4

Version : 0112

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Voltmeter sensor

1 Généralités

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1 General

This sensor measures instantaneous voltage values up to ± 24 V.
The measurements are made by a differential input with safety banana jacks.

2 Use

This sensor operates with the Tooxy console in standalone mode (without computer), or in connected mode (with computer) after being connected to the Tooxy console (4mm diameter bushings).

2.1 Connection to the Tooxy console

The sensor is connected using the connector on the back of the housing:

- Bring the sensor in line with the selected measurement channel
- Insert the sensor along the console guide rail
- follow the Tooxy console guide system
- Push all the way in

2.2 Type of measurement

The sensor measures the instantaneous value of the voltage at its terminals.

2.3 Choice of the range

The two available ranges are ± 12 V and ± 24 V.
A button on the sensor housing is used to modify the sensor range.

To modify the sensor range

2.4 Operation in standalone mode

a. Start up

- connect the sensor
- switch the console on

b. Display

An instantaneous value of the voltage is displayed directly on the display screen on the outside of the housing.

2.5 Operation in connected mode

a. Start up

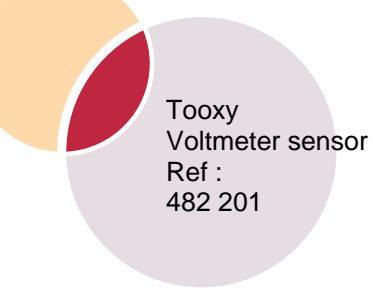
- connect the sensor to the console
- switch the console on
- connect the console to the computer

b. Display

Measurements are displayed directly on the computer display screen when an acquisition is made or using the "display" tab.

3 Technical characteristics

Measurement range:	- range ± 12 V	from -12 V to $+12$ V
	- range ± 24 V	from -24 V to $+24$ V
Resolution	- range ± 12 V	6 mV
	- range ± 24 V.	12 mV
Precision	- range ± 12 V	$\pm (1\% + 2\text{LSB})$ of the value
	- range ± 24 V.	$\pm (1\% + 2\text{LSB})$ of the value
Passband	≥ 10 kHz	
Protection	140 VAC / ± 200 VDC	

Tooxy
Voltmeter sensor
Ref :
482 201

4 After-Sales Service

This material is under a two year warranty and should be returned to our stores in the event of any defects.

For any repairs, adjustments or spare parts, please contact:

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SA capital 3 233 762 € - Siren R.C.S. B 387 901 044 - Siret 387 901 04400017

Direct connection for technical support

A team of experts at your disposal from Monday to Friday (opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediately to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

Free service *

+ 33 (0)2 32 29 40 50**

* Call cost not included

** Only for call from foreign countries