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Faune terrestre

## *Environment*

*Fauna*

Ref :  
525 001

Français – p 1

English – p 3

Version : 6010

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### *Insect aspirator*

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## 1 Introduction - General description

This insect collector is specially designed for a safe use by each student.

Two washable and interchangeable mouthpieces are supplied for preventing any risk of contamination. A mechanical filter, essential during use, prevents the passage of insects into the suction tube.

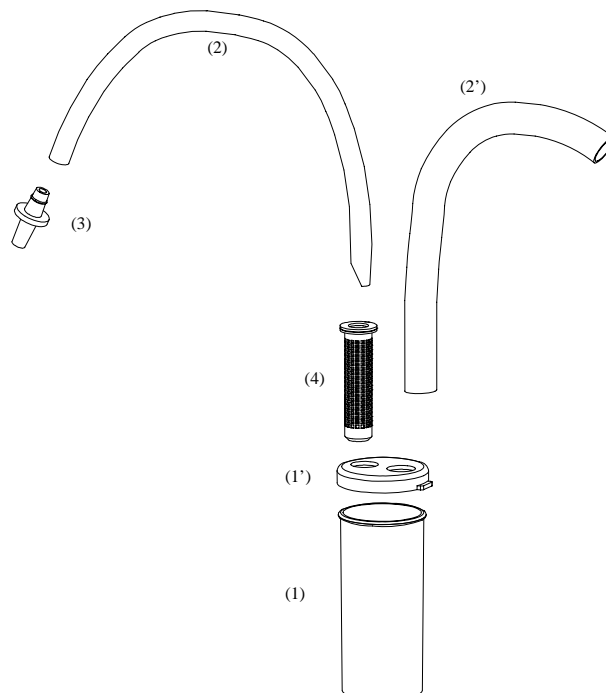
It allows capturing small insects without damaging them. It is suitable for capturing and sorting the **drosophila** to be selected for your experiments in genetics.

It can be fully dismantled and washed for cleaning and disinfection.

## 2 Composition

- A container (Ø35 mm x 80 mm) **(1)**
- A cover with hole **(1')**
- Two tubes: one for sucking and the other for placing near the insects to be captured **(2)** and **(2')**
- Two washable mouthpieces for preventing any risk of contamination between the users **(3)**
- A mechanical filter, non removable for preventing any insects from being sucked in accidentally **(4)**.

## 3 Set-up

## 4 Procedure

Mount the collector as shown above.

Place the larger diameter tube near the insect to be captured. Suck via the mouthpiece; when the insect is sucked, it remains captive in the collector container.

## 5 Suggestions for usage

- Capturing small living insects without damaging them.
- Sorting the anesthetised **drosophila** to be selected for crossing the strains.

## 6 Storage

Dismantle the collector and clean the different parts with disinfectant. Allow to dry before storage.

## 7 After-Sales Service

This material is under a two year warranty and should be returned to our stores in the event of any defects.

For any repairs, adjustments or spare parts, please contact:

**JEULIN - TECHNICAL SUPPORT**  
**Rue Jacques Monod**  
**BP 1900**  
**27 019 EVREUX CEDEX FRANCE**  
**+33 (0)2 32 29 40 50**

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SA capital 3 233 762 € - Siren R.C.S. B 387 901 044 - Siret 387 901 04400017

## Direct connection for technical support

A team of experts at your disposal from Monday to Friday (opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediatly to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

**Free service \* :**  
**+ 33 (0)2 32 29 40 50**

\* Call cost not included