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Measurement and instrumentation

Ref :
291 194

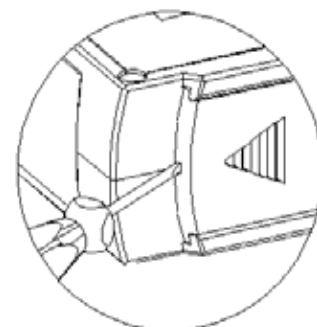
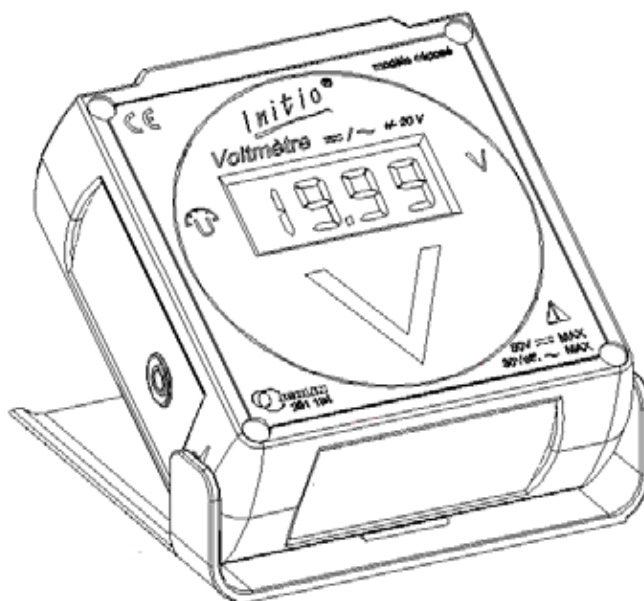
Français – p 1

English – p 3

Version : 6010

Voltmètre A/C Initio[®]

Voltmeter A/C Initio[®]

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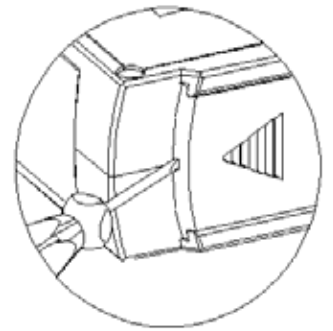
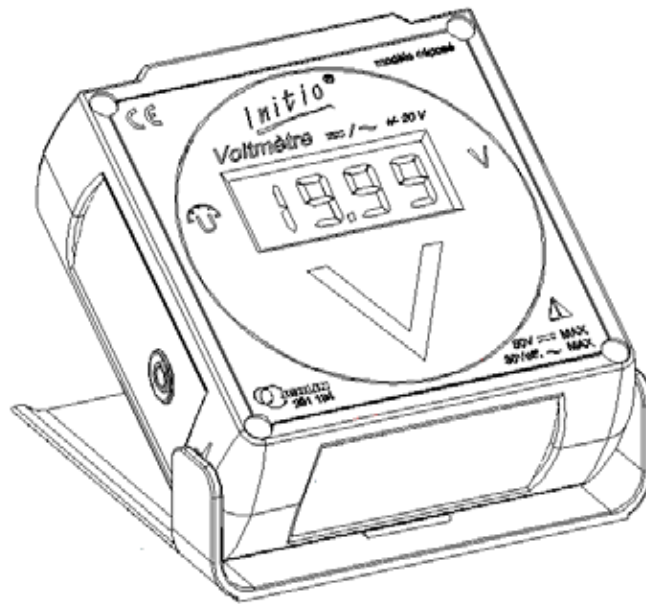
1 Technical characteristics

- Measuring range :
 - +/- (0.00 to 19.99) V continuous
 - (0.070 to 19.99) V effective
- Resolution :
 - 10 mV
- Precision :
 - +/- (0.5% VL + 1 UR) continuous
 - +/- (1.5% VL + 4 UR) alternative
- Voltage :
 - 50 V continuous and 30 V effective
- Protection :
 - up to 230 V effective
- Power supply :
 - 1 battery, 9 V 6F22 type

2 Functioning

1 – Set up of the device is carried out by connecting a Ø4mm banana lead to the red socket.

2 – The \sim/\neq selection is automatic.

- if the effective value of the signal goes over 700 mV, the device will display an effective voltage value (alternative mode).
- for an effective value less than 700 mV, the device will display a mean voltage value (continuous mode).

3 Change of battery

When the aide 'Bat' appears in the top left of the display, it is recommended to change the battery. In order to open the battery casing, use a small screwdriver, inserting it into the notch on the edge of the door, and lifting it as indicated in the drawing on the first page.

4 Resetting

Although stable in the workshop, it is possible to return the value to ZERO with the help of a potentiometer situated at the base of the battery unit. For this, insert a small screwdriver, in the hole on the right, and turn until 0.000 is obtained by having previously lined up the two inputs between them. This adjustment corresponds to the offset adjustment, and will displace the measures, more or less, to the same value.

5 After-sales service

The device is under a 2-year guarantee, it must be sent back to our workshops.

For any repairs, adjustments or spare parts please contact:

JEULIN – TECHNICAL SUPPORT
Rue Jacques Monod
BP 1900
27 019 EVREUX CEDEX FRANCE
+33 (0)2 32 29 40 50

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SA capital 3 233 762 € - Siren R.C.S. B 387 901 044 - Siret 387 901 04400017

Direct connection for technical support

A team of experts at your disposal from Monday to Friday (opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediately to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

Free service * :
+ 33 (0)2 32 29 40 50

** Call cost not included*