Son

Sound

**Ref :
302 136**

Français – p 1

English – p 3

Version : 6010

Ampli-sons

Ampli-sons

1 Objectifs

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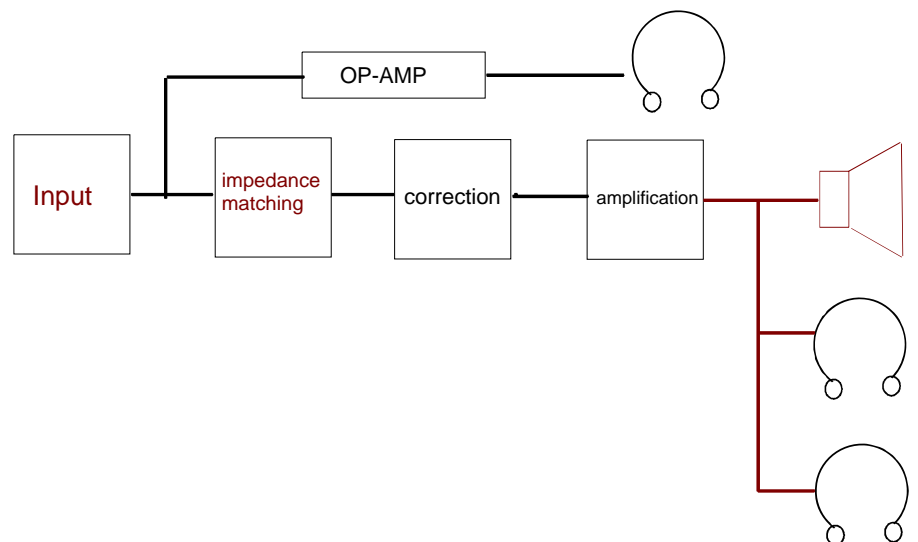
1 Objectives

- Amplification by means of a single op-amp wired as a non-inverting amplifier
- Power amplification with:
 - impedance matching
 - treble and bass correction
- Solving connection problems between the audio instrument and the measuring instrument

2 Principle

The input signal can be fed via a male jack input or via two male banana plugs. The signal is then amplified in two parallel-mounted independent stages having specific outputs:

- an op-amp wired as a non-inverting amplifier having a gain of 10–100 and fitted with a high-impedance headphones output
- a power amplifier (10W) with one loudspeaker output and two high-impedance headphones outputs.

3 Technical data

- Dimensions : 160 x 90 x 120 mm
- Mass : approx. 0.3 kg
- Power supply : -15V; 15V
- Protection : against voltage surges or polarity inversion
- Output voltage : - op-amp: +15V;-15V
- power amplifier: 0-10V
- Output power : - power stage: 10W
- op-amp: a few mW ; gain = 10 to 100
- Correction : - 3 Hz < treble < 16 kHz
- Bass < 400 Hz
- Bandwidth : 40 Hz to 16 kHz
- Input connector : - female jack (dia. 6.35 mm)
- banana plugs (dia. 4 mm)
- Output connectors : - op-amp: one female stereo jack (dia. 3.5 mm)
- power amplifier : banana plugs (dia. 4 mm)
- 2 female stereo jacks (dia. 3.5 mm)

4 After-Sales Service

This material is under a two year warranty and should be returned to our stores in the event of any defects.

For any repairs, adjustments or spare parts, please contact:

JEULIN - TECHNICAL SUPPORT
Rue Jacques Monod
BP 1900
27 019 EVREUX CEDEX FRANCE
+33 (0)2 32 29 40 50

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SA capital 3 233 762 € - Siren R.C.S. B 387 901 044 - Siret 387 901 04400017

Direct connection for technical support

A team of experts at your disposal from Monday to Friday (opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediatly to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

Free service * :
+ 33 (0)2 32 29 40 50

* Call cost not included